SPECTRO-UV[®]

**QZ-Series
UV-A Magnifier Lamps
OPERATOR'S MANUAL**

SPECTRO-UV[®]

www.Spectro-UV.com
4 Dubon Ct., Farmingdale, NY 11735
866-230-7305

PRINTED IN U.S.A.

AM11002ML-3
ISSUED: OCTOBER 2023

SPECTRO-UV® QZ-Series
UV-A Magnifier Lamps

 **DANGER**

Ultraviolet radiation emitted from this product. Avoid exposure. **ALWAYS WEAR PROTECTIVE CLOTHING. EXPOSURE MAY CAUSE PREMATURE AGING OF THE SKIN AND CANCER. ALWAYS WEAR PROTECTIVE EYEWEAR; FAILURE TO DO SO MAY RESULT IN SEVERE BURNS OR LONG TERM INJURY TO EYE.** Never look directly into the lamp. Exposure can cause eye and skin allergy and allergic reactions. Medications or cosmetics may increase your sensitivity to ultraviolet radiation. Consult physician before operating this product if you are using medications or have a history of skin problems or believe yourself especially sensitive to sunlight.

FOR LABORATORY USE ONLY

UNPACKING AND INSPECTION

1. Carefully unpack and inspect the lamp for possible damage in shipment. If any damage is noted, immediately notify the carrier and supplier before attempting to use the lamp.
2. Save the shipping carton and packing materials for future storing and shipping of the lamp.
3. Check the lamp's electrical performance as soon as possible according to the GENERAL FUNCTIONAL CHECKS outlined below.

 **WARNING**

Please read the instructions for this lamp carefully before using it. **Never use this lamp in any manner not specified in these instructions. Do not use if a tube, cord or plug is not fully serviceable. We recommend using Spectro-UV® UV-absorbing protective eye and face wear. See Replacement Parts and Accessories for model numbers. This eye and face wear also improves contrast between the fluorescent area and the background and eliminates "blue haze" interference.**

GENERAL FUNCTIONAL CHECKS

1. Connect the plug to a power outlet that complies with the electrical requirements specified on the lamp's label.
2. The power supply cord is the disconnect for this unit.
3. Press the switch to the "I" position to operate the UV-A long wave tubes. Press the switch to the "II" position to operate the white lights. The "O" position turns the lamp off.
4. The tubes should become energized and emit a steady glow of light. The light may initially flicker, especially if the lamp is cold, but should stabilize after a few seconds.

OPERATOR SERVICING

The following servicing instructions are intended for limited maintenance of the lamp. Do not perform any servicing other than that contained in these instructions. Return the lamp to the factory for other maintenance not covered in this section.

WARNING
To avoid the risk of electrical shock, always unplug the lamp from its power source before cleaning or servicing.

HOUSING AND MAGNIFIER LENS MAINTENANCE

The polycarbonate lamp housing protects the internal parts. Immediately clean spills from the housing and magnifier lens with water and a mild detergent and wipe dry with a soft cloth. Do not allow any liquids to leak into electrical components. Periodically clean the lens with a glass cleaner and a soft cloth.

TUBE AND REFLECTOR MAINTENANCE

To clean the tubes and/or reflector, see the "TUBE REPLACEMENT" section. Using a soft cloth, the reflector may be cleaned with a glass cleaner, while the tubes may be cleaned with a mild detergent solution. Wipe dry with a soft cloth.

TUBE REPLACEMENT

When replacing a tube, unplug the lamp from its power source before proceeding.

1. Lay the lamp housing on a flat, level surface with the tubes facing up.
2. Gently rotate each tube (one at a time) until the metal contacts at each end line up with the small rectangular openings. The tube may now be easily removed from its sockets.
3. Install the new tube by first aligning the contacts at the end of each tube with the small rectangular openings at each end of the sockets.
4. Turn one tube at a time gently to rotate about 90° for the tube to be locked into place by the "Safety Slots." *Safety Slots* ensure proper contact by aligning each tube, provide insulation and protects from electrical shock.
5. If installation of the new tube fails to yield correct lamp operation, check to make sure the two pins at each end of the tube are properly connected with the two copper contacts in each socket.

NOTE: The tubes in all QZ-Series lamps are wired in series. Therefore, if one tube fails the other tube of the same wavelength will not light, even though it may be in good operating condition.

ENVIRONMENTAL SPECIFICATIONS

The Spectro-UV QZ-Series lamps are designed to be safe under the following conditions:

- Indoor use;
- Altitude up to 2,000 m (6,562 ft.);
- Temperature 5°C to 40°C (41°F to 104°F);
- Maximum relative humidity 80% for temperatures up to 31°C (88°F), decreasing linearly to 50% relative humidity at 40°C (104°F);
- Mains supply voltage fluctuations not to exceed $\pm 10\%$ of the nominal voltage;
- Installation Category II;
- Pollution Degree 2.

ELECTRICAL SPECIFICATIONS

MODEL	WAVELENGTH	NO. TUBES	AMPS
QZ-22	UV-A/WHITE	2/2	0.2
QZ-22B	UV-A	4	0.4

All Domestic models are 120V/60Hz.

All European (/F) and British (/FB) models are 230V/50Hz. CE compliant.

All Australian (/FA) models are 240V/50Hz.

All Japanese (/J) models are 100V/50-60Hz.

Housing dimensions W x L x H: 5.75 in x 7 in x 2.25 in (14.6 x 17.8 x 5.7 cm).


Tubes are 4 Watt. The UV-A tubes are integrally filtered.

Long Wave (UV-A): 365nm

Front View

Back View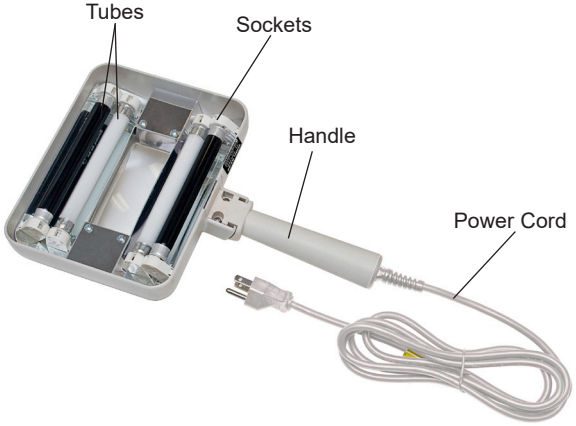

REPLACEMENT PARTS AND ACCESSORIES

Description

Tube, Long Wave UV-A
Tube, White Light
Spectacles, UV-Absorbing
Goggles, UV-Absorbing
Face Shield, UV-Absorbing

Part No.

BLE-220B
BLE-95D
UVS-30
UVG-50
UVF-80

WARRANTY

The warranty policy for the QZ-Series lamps is provided on the Certificate of Limited Warranty enclosed separately with each unit.

NOTE: For assistance of any kind, including help with a lamp under warranty, contact the Customer Service Department at Spectro-UV. In the U.S. and Canada, call toll-free 1-866-230-7305. Include the model and serial numbers of the unit and the date of purchase. Shipping instructions will be provided if necessary.

If an estimate of charges for nonwarranty work or other service work is required, a quote will be furnished upon evaluation of the unit. Service work will not be performed without customer approval.

SHIPPING

Carefully pack the lamp in the original shipping container and packing materials. Insure it for full value and ship it prepaid to the factory.

LAMP CONTAINS MERCURY
Manage in Accord with Disposal Laws
See: www.lamprecycle.org

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